



# Mass Spectrometry Solutions

WWW.MKSINSTITUTE.COM

## Cirrus™ 3

### POWERFUL AND VERSATILE ATMOSPHERIC GAS ANALYSIS

Designed for researchers and engineers who require fast, sensitive, reliable monitoring of gases at atmospheric pressure, the Cirrus™ 3 system provides superior analytical results in a compact, powerful, versatile platform for the lab or production facility.

Building on proven Cirrus Quadrupole Mass Spectrometry technology, the Cirrus 3 combines state-of-the-art performance in compact benchtop and rack-mounted formats. With a heated capillary inlet and fast scan speed (250 points per second), gas compositions can be tracked over wide dynamic range (ppb to percentage levels) more reliably.

Enhancements in capillary and oven design ensure precise temperature control and stability, allowing you to obtain high quality data with greater confidence in your results. The versatile Cirrus 3 system with its automatable software control can be optimized and customized for a variety of applications, from industrial processes to research applications.

Featuring automatic start-up and shut-down routines and built-in vacuum and heater interlocking for system protection, Cirrus systems are easy to install and operate. Cirrus 3 systems are manufactured from quality, field-proven materials to ensure reliability and maximize up-time. Particular consideration has been paid to ease of service and maintenance.

### Features & Benefits

#### Reliable Performance

- Monitors multiple gas species over a wide dynamic range from trace levels (ppb) to high abundance (percentage levels)
- Fast response – 250 data points per second, and capillary inlet heated up to 300°C
- Precise and stable user-defined temperature control with enhanced oven and capillary designs
- Detect gases up to 300 amu. High resolution options (1-100HP)

#### Versatile and Easy to Use

- Recipe-driven Process Eye™ Professional supports gas monitoring and quantitation applications
- Customizable 'Recipe-based' software provides automated operation and calibration

- Fully network compatible via direct Ethernet interface
- Compact benchtop and rack-mount formats ideal for laboratory and manufacturing environments
- Modular design for ease of maintenance and service
- Multiple inlet options for sampling different gas conditions:
  - Fused silica (up to 150°C) and stainless steel (up to 300°C) capillaries
  - Low flow capillaries
  - Multi-stream inlet (4-, 8-, and 16-way)
  - Automated variable pressure inlet control (option)



## Applications

- Monitoring trace contaminants in process gases
- Gas purity and manufacturing QC
- Catalyst studies – Chemisorption
- Thermal analysis – TGA, DTA, DSC
- Fuel cell monitoring and development
- Fermentation and exhaust monitoring
- Heat treatment/furnace monitoring
- Membrane studies
- Glove box gas monitoring
- Lamp manufacturing
- Freon detection and identification

## Introduction

Cirrus systems are based on Quadrupole Mass Spectrometry, widely acknowledged as a preferred platform technology for many atmospheric pressure gas analysis applications. The Cirrus 3 system, with its improved temperature control and compact design, provides high levels of performance and robustness in space-efficient and flexible configurations that are ideal for a wide range of analytical applications.

Precision engineered to provide fast and reproducible data, Cirrus 3 can accurately detect and quantify gases and gas mixtures, from one or more sampling locations with a single instrument. Gas composition can be monitored over a wide dynamic range, from percentage levels down to the sub ppm range.

Cirrus 3 is an ideal choice for atmospheric gas analysis whether you work in a research setting or a manufacturing environment. By combining Process Eye Professional software with MKS technical support, Cirrus can be integrated, optimized, and automated to your precise specifications. Ultimately, users across your organization realize greater benefits of the power of quadrupole mass spectrometry.

## Quadrupole Analyzer

At the heart of every Cirrus 3 system is a precision-built quadrupole analyzer incorporating a closed ion source, a triple mass filter, and a dual (Faraday and Secondary Electron Multiplier) detector system. This analyzer configuration is selected to optimize sensitivity and long-term performance stability.

## Heating and Vacuum System

The Cirrus 3 features a number of advancements in design that aid stability and reproducibility of results. The vacuum chamber (with quadrupole analyzer) and inlet interface assembly is housed within a precision engineered stainless steel oven, providing low heat sink characteristics. This offers high temperature uniformity and eliminates potential cold spots that could otherwise cause undesirable sample condensation and variance in analyses. The oven design enables user-defined temperature control of the oven and capillary, allowing for finer analytical control.

The quadrupole analyzer itself is contained within a stainless steel vacuum chamber which is pumped by a high compression turbo molecular/diaphragm pump combination. The vacuum system is backed by an internal 4-stage diaphragm pump to maintain the integrity of the compact design. The entire Cirrus 3 vacuum chamber can also be baked to reduce residual gas background species and to minimize any memory effects.

## Gas Inlet

An essential feature of any gas analyzer is the ability to not contaminate or alter the gas sample in any way. The Cirrus 3 inlet assembly incorporates either fused silica or stainless steel capillaries that can be heated to a stable user-defined temperature. The inlet assembly allows for easy replacement of capillaries, while the optimized insulation maximizes stability and reduces heat radiation.

The low volume and surface area of the assembly serves to maximize response speed while minimizing memory effects. Cirrus 3 systems can also be configured with inlets for multi-stream sampling, stainless steel capillaries for resistance to fluorine-based compounds, and a pressure controller inlet (incorporating a gas independent capacitance manometer) to allow sampling from sources that vary from the nominal 1 bar inlet requirement. The Cirrus 3 vacuum system utilizes a high compression turbo molecular pump, so light gases such as hydrogen and helium can be sampled with no additional expensive pumping requirements.

## Maintenance and Access

The Cirrus 3 internal oven has a removable cover allowing easy access to the inlet interface, vacuum chamber, and analyzer ion source. This is particularly helpful for routine maintenance, such as filament and capillary replacement. A hot cathode (ion) gauge is incorporated for independent vacuum pressure measurement and to provide an interlock signal for protection of the mass spectrometer. A temperature sensor also ensures that the electron multiplier detector cannot be switched on at high temperatures, preserving its life. The Cirrus 3 is designed with a lubricant-free pumping system. No elastomer seals are used in the sample inlet system or in the high vacuum region of the system.



## Process Eye™ Professional Software

Cirrus 3 is operated using proven Process Eye Professional software, a recipe-driven platform that communicates with the system over a TCP/IP network. Process Eye Professional is designed for use with the latest Microsoft® operating systems including 32 bit or 64 bit Windows® 7 and 10.

Features and Benefits include:

- Data presented in units relevant to the application
- Fully automated operation and calibration
- New workflow "Recipes" for quantitative applications
- User-configurable alarms and warnings
- Can be configured to track data from other process sensors (temperature, pressure, flow, etc.)

## Options

- Multiple gas stream inlet (4-, 8- or 16-way)
- Corrosive gas sampling version
- Regulated ion source pressure option to compensate for variation in inlet pressure and gas composition
- High mass resolution (1-100 amu HP)



Cirrus 3 rack-mount configuration

## Specifications

### Dimensions & Weight

Benchtop System  
Rack-mount System

593mm L x 291mm W x 470mm H, 35 Kg  
577mm L x 482mm W x 266mm (6U) H, 35 Kg

### Mass Range

1-100 amu; 100HP, 1-200 amu and 1-300 amu options

### Detection Limits

Gas dependent, typically < 100 ppb for non-interfering species

### Electron Energy & Emission Current

Operator variable

### Operating Temperature and Humidity Range

5°C to 35°C, 20% to 80% RH (non-condensing)

### Oven Temperature

180°C for bakeout. User-definable for operation – typically 80°C

### Capillary Inlet

2.0m long with ¼" Swagelok® end connection, heated to 150°C (optional heating to 300°C, with stainless steel only). Standard fused silica and optional stainless steel. User-definable temperature control.

### Gas Consumption

20 ml/min (Lower uptake rate capillaries options are available)

### Sample Pressure

1 bar nominal

### Pumping System

High compression turbo-molecular pump with internal 4-stage diaphragm backing pump standard. Corrosive gas tolerant pumping with all internal pumps optional.

### Vacuum Hardware

60 l/s turbo molecular pump with high conductance analyzer housing, inlet system, automated vacuum controller (RVC) completely interlocked and integrated.

### Automated Inlet Pressure Controller

A 640 (Baratron® capacitance manometer based) automated inlet pressure controller is available as an option

### Recommended PC Spec

Microsoft 32 bit or 64 bit Windows 7 or 10

### Computer Interface

1 x RJ45 Ethernet LAN port required

### I/O Capability

4 analog inputs (-11 to +11 volt, 22 bit)  
2 analog outputs (0-10 volt, 12 bit)  
16 TTL digital I/O

### Power

Universal mains input 100 – 240 VAC / 50 – 60 Hz  
Power required: 700W maximum (during bakeout); 300W typical

### Compliance

CE



# Ordering Information

## Rack-mount Instrument Product Ordering Codes:

### Non-process

Cirrus 3 (100 amu, non-process): 473-12E-A30  
 Cirrus 3 (200 amu, non-process): 473-22E-A30  
 Cirrus 3 (300 amu, non-process): 473-32E-A30

### Process\*

Cirrus 3 (100 amu, process): 473P-12E-A30  
 Cirrus 3 (200 amu, process): 473P-22E-A30  
 Cirrus 3 (300 amu, process): 473P-32E-A30

## Benchtop Instrument Product Ordering Codes:

### Non-process


Cirrus 3 (100 amu, non-process): 474-12E-A30  
 Cirrus 3 (200 amu, non-process): 474-22E-A30  
 Cirrus 3 (300 amu, non-process): 474-32E-A30

### Process\*

Cirrus 3 (100 amu, process): 474P-12E-A30  
 Cirrus 3 (200 amu, process): 474P-22E-A30  
 Cirrus 3 (300 amu, process): 474P-32E-A30

\* Process packages include on-site applications start-up assistance.

For other options such as filament type, corrosion resistance, inlet pressure control, lower inlet flow capillaries and capillary material and temperature, contact a local MKS sales office for applications assistance.

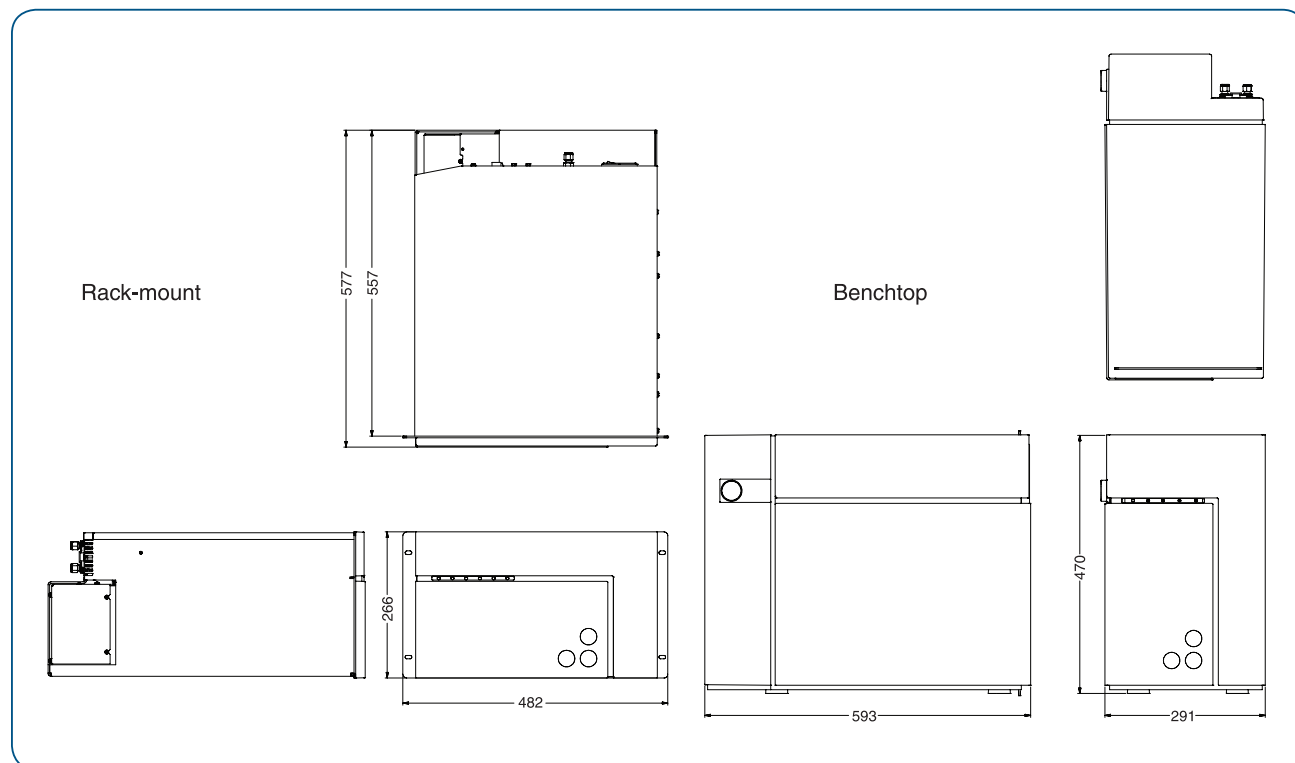

## Dimensional Drawing —

Note: Unless otherwise specified, dimensions are nominal values in millimeters.



### MKS Instruments, Inc. Global Headquarters

2 Tech Drive, Suite 201  
 Andover, MA 01810  
 Tel: 978.645.5500  
 Tel: 800.227.8766 (in USA)  
 Web: www.mksinst.com

### MKS Instruments, Inc. Mass Spectrometry Solutions

3635 Peterson Way  
 Santa Clara, CA 95054  
 Tel: 408.750.0300

### MKS Instruments, Inc. Mass Spectrometry Solutions

Cowley Way  
 Crewe, Cheshire CW1 6AG  
 Tel: +44.1270.250150